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ENGLISH

A. Parts and Components

The following illustrations show the names of each component, button and switch on the microscope.

1. Optical switch
2. Power indicator
3. Power & Zoom in & Menu
4. LCD screen
5. Function (Right)
6. Playback (Left)
7. USB port
8. SD card slot
9. Strap eyelet
10. Shutter / OK
11. TV-out (optional)
12. Battery compartment cover
13. View case (optional)
14. Lens
15. Tripod / screw socket
16. LED
17. Magnification switch

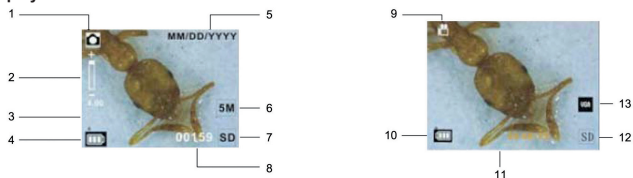
B. Quick start

- Installing microscope batteries (not included):
 1. Slide open the battery compartment cover.
 2. Insert the batteries. Be sure to load the batteries with + and – terminals as shown on the microscope.
 3. Close the battery compartment (slide it up).
- Inserting a SD card (not supplied):
 1. Turn off the microscope.
 2. Insert the SD card as shown. Push the card into the SD card slot.
 3. To remove the card, push it and it will release.
- Turning on the microscope:
 1. Press the POWER button as shown in the illustration.
 2. The indicator will light up and the LCD screen will turn on.
 3. Press the POWER button again to shutdown the microscope.

C. Display information

- | | |
|---------------------------|---|
| 1. Capture mode indicator | 8. Memory capacity (remaining number of storable pictures) |
| 2. Zoom indicator | 9. Video mode indicator |
| 3. Single picture icon | 10. Battery capacity |
| 4. Battery capacity | 11. Memory capacity (remaining record time of storable video) |
| 5. Date | 12. SD card |
| 6. Picture size | 13. Video size |
| 7. SD card indicator | |

D. Using the microscope in PC cam mode

Note: Please install the microscope driver before you use the microscope in PC cam mode.

Setting up PC cam

- Turn the microscope power on and press the FUNCTION button to select the setup function. Press the OK button to confirm. Press the RIGHT button to select USB connection. Select the PC cam. Press the OK button to confirm.
- Connect the microscope via the USB cable to a computer.
- There are many methods to view your image in PC cam mode if you have video software. For example: Double-click "My Computer", and then click "Standard Camera". You can live view the PC cam mode now.

For more functions check the full manual on CD.

In this manual you can find the specifications and advanced functions of the device.

Safety precautions:

To reduce risk of electric shock, this product should ONLY be opened by an authorized technician when service is required. Disconnect the product from mains and other equipment if a problem should occur. Do not expose the product to water or moisture.

Maintenance:

Clean only with a dry cloth. Do not use cleaning solvents or abrasives.

Warranty:

No guarantee or liability can be accepted for any changes and modifications of the product or damage caused due to incorrect use of this product.

General:

Designs and specifications are subject to change without notice.

All logos brands and product names are trademarks or registered trademarks of their respective holders and are hereby recognized as such.

Keep this manual for future reference.

Attention:

This product is marked with this symbol. It means that used electrical and electronic products should not be mixed with general household waste. There is a separate collections system for these products.

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Copyright ©
NEDIS B.V.