MANUAL (p. 2)
USB Microscope

ANLEITUNG (S. 4)
USB-Mikroskop

MODE D'EMPLOI (p. 6)
Microscope USB

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
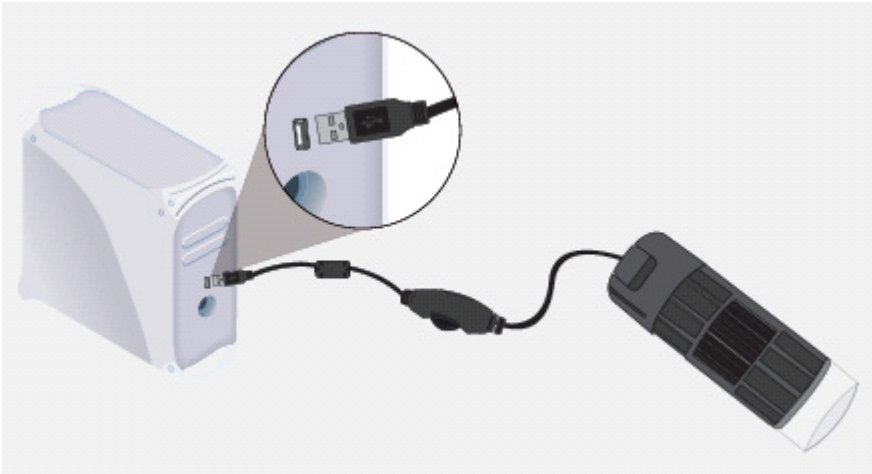
ENGLISH

1. Connecting the microscope

- Plug the USB cable into the computer. The device will be automatically detected in Windows® XP or later versions.

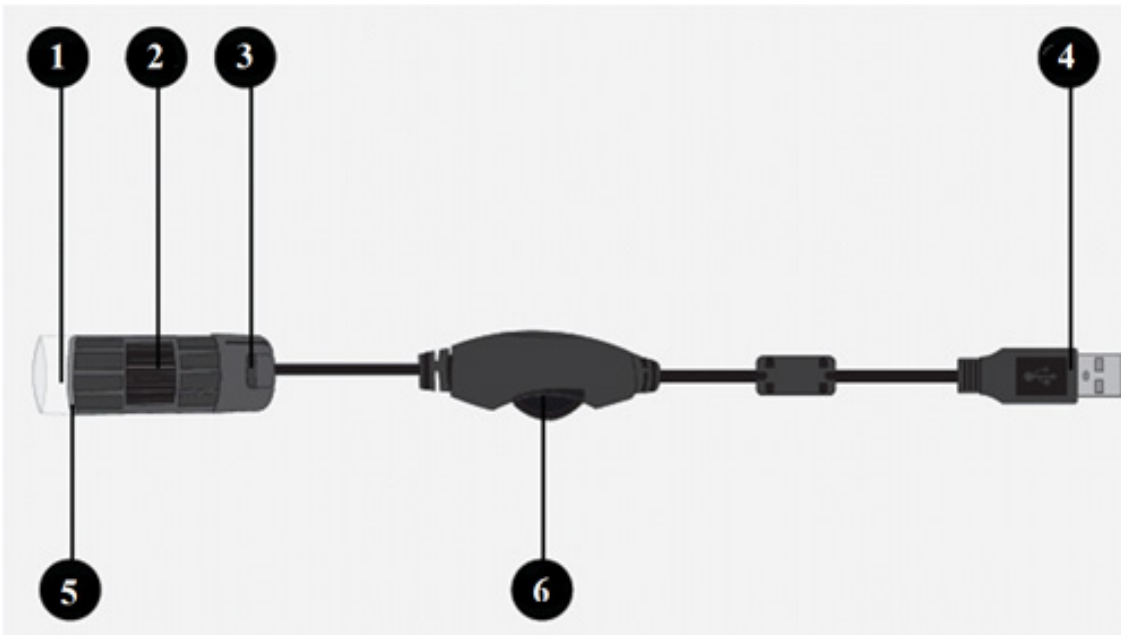- The device is compatible with any editing software.

- | | |
|-------------------|--|
| 1. Lens: | Place over the desired object to watch |
| 2. Focus: | Turn left or right to set focus |
| 3. Snapshot: | Press to get a snapshot of the current image (does not work with some software versions) |
| 4. USB connector: | Used to connect the microscope to the PC |
| 5. LED lights: | Turn on when needed |
| 6. Light control: | Move left or right to dim or turn the light on/off |

3. Software installation

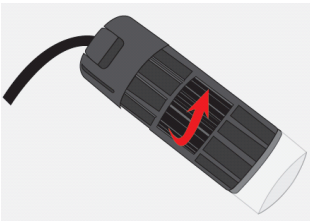
- The microscope has some viewing software, however you can also use other webcam software with webcam functions (such as MSN messenger, Skype etcetera).
- To install the software, copy only the “amcap2a” software file to your desktop.

4. How to use the microscope


- After connecting the microscope to the computer, run the “amcap2a” software.
- A new window will appear showing the microscope image.
- Place the microscope over the desired viewing area.

- Set the focus.

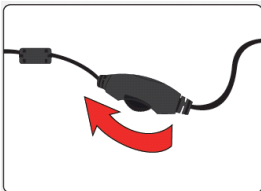- In case you need additional light, use the light control to turn the LED's on.

- When the microscope is not in use, it is advisable to unplug it and re-connect it the next time.
- ⚠ If the image has not been displayed and another device is also in use (such as a webcam), you may need to select the “Devices” menu and change the input to USB 2.0 camera.

Safety precautions:

CAUTION
RISK OF ELECTRIC SHOCK
DO NOT OPEN

To reduce risk of electric shock, this product should **ONLY** be opened by an authorized technician when service is required. Disconnect the product from mains and other equipment if a problem should occur. Do not expose the product to water or moisture.

Maintenance:

Clean only with a dry cloth. Do not use cleaning solvents or abrasives.

Warranty:

No guarantee or liability can be accepted for any changes and modifications of the product or damage caused due to incorrect use of this product.

General:

Designs and specifications are subject to change without notice.

All logos brands and product names are trademarks or registered trademarks of their respective holders and are hereby recognized as such.

Keep this manual and packaging for future reference.

Attention:

This product is marked with this symbol. It means that used electrical and electronic products should not be mixed with general household waste. There is a separate collections system for these products.

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Copyright ©
NEDIS B.V.